



---

## **NYX1-512 Datasheet**

512 Electrode Neural Recording ASIC

Product Datasheet (Version 1.4)

Science Corporation  
300 Wind River Way, Alameda, CA 94501


[www.science.xyz/technologies/ecosystem](http://www.science.xyz/technologies/ecosystem)  
© Science Corporation

# Table of Contents

<a href="#">Table of Contents</a>	1
<a href="#">Summary</a>	2
<a href="#">Key Features</a>	2
<a href="#">Applications</a>	2
<a href="#">Architecture</a>	3
<a href="#">Summary Table</a>	4
<a href="#">Electrical Interface</a>	6
<a href="#">Digital I/Os</a>	7
<a href="#">Command Structure</a>	9
<a href="#">The Nixel Array</a>	10
<a href="#">Electrode Input Arrangement</a>	11
<a href="#">Recording Parameters</a>	12
<a href="#">Troubleshooting</a>	12
<a href="#">Recording Configuration</a>	13
<a href="#">Recording Setup Sequence</a>	13
<a href="#">Reset</a>	13
<a href="#">Enable Channels</a>	13
<a href="#">Configure LNA Inputs</a>	13
<a href="#">Set Sample Rate and Bit Depth</a>	14
<a href="#">Configure Reference Routing</a>	17
<a href="#">Initiate Start Sequence</a>	17
<a href="#">Stop Recording</a>	17
<a href="#">Pin Specifications</a>	18
<a href="#">On-Chip Registers</a>	23
<a href="#">AnaDrive</a>	23
<a href="#">Nixel</a>	25
<a href="#">ADC</a>	26
<a href="#">DigCont</a>	27
<a href="#">Unused</a>	30
<a href="#">Packaging</a>	31
<a href="#">Contact Information</a>	33

## Summary

The NYX1-512™ chip is a low-noise and low-power mixed-signal programmable neural recording ASIC with 512 electrodes and 16-bit data path. It has an array of 256 differential record channels with integrated low-noise amplifiers (LNAs) and analog-to-digital converters (ADCs) that can record 512 electrodes in parallel at sampling rates up to 32 kHz with ADC resolutions up to 12-bit. The chip has a uniform array of 256 fully differential identical active neural recording elements, called nixels, acting as electrical pixels for neural recording applications. The NYX1-512 chip has a tiny chip-scale package measuring 4 mm x 4 mm, suitable for ultra compact head stages and high-density multi electrode-array (MEA) neural recording applications.

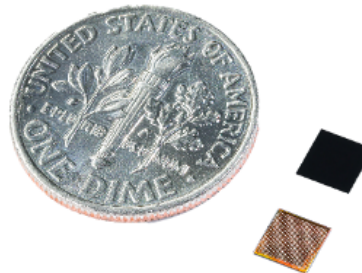

## Key Features


- 512 recording electrodes in a uniform array.
- 256 fully differential neural recording elements (nixels).
- User-selectable analog polarity for input and reference electrode selections.
- AC coupled LNA with in-nixel digitization with digital readout.
- Programmable nixel gain and bandwidth (BW) selection for low noise.
- Input referred noise below 5  $\mu\text{V}$  rms in the high gain mode.
- High- and low-gain modes for spike and local field potentials (LFP) recording.
- User-selectable reference generation and distribution.
- Simultaneous spike and LFP recording up to 32 kHz sampling rate.
- Programmable ADC resolution from 10 to 16-bit.
- Impedance measurement capability for all 512 electrodes.
- Low-power and low-voltage design (1.2 V digital and 2.5 V analog).
- Wafer level chip-scale package (WLCSP): 4 mm x 4 mm (Cu bumps).

## Applications

- Chip-scale ultra-compact neural recording solutions for high-density MEAs.
- Ultra-compact high-density head stages with digital control and readout.
- Active MEAs for “Smart-Dish” type in vitro neural recording applications.
- Chip-scale MEAs for “Neurons-on-Chip” type neural recording applications.
- Low-noise multi-channel scientific data acquisition systems.

## Architecture

The NYX1-512 chip is composed of the nixel array, analog drive, and digital drive.

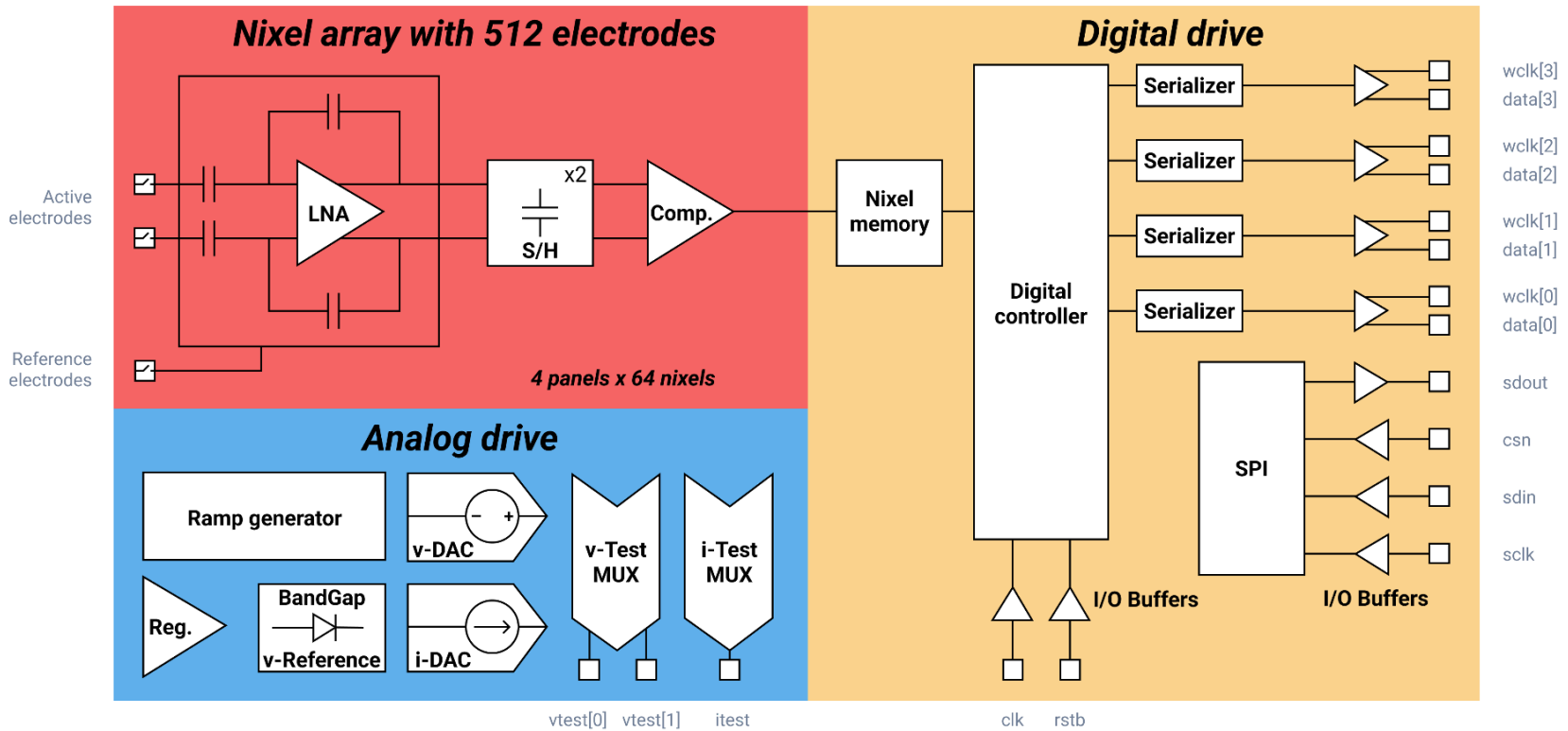

Summary Table


Power supplies and returns (grounds)	Analog power		vdda_bg	Bias generator supply	2.5 V
			vdda_lna	Nixel LNA supply	
			vdda_cmp	Nixel comparator supply	
	Analog ground		vssa_bg	Bias generator ground	0.0 V
			vssa_lna	Nixel LNA ground	
			vssa_cmp	Nixel comparator ground	
			sub	Substrate, ground	
Digital power		dvdd2p5	Level shifter supply	2.5 V	
		dvdd	Core and I/O supply	1.2 V	
Digital ground		dvss	All digital block, ground	0.0 V	
Power dissipation	32 KHz sampling 12-bit ADC resolution 160 MHz clk	Analog	≤ 22mW		
		Digital	≤ 8mW		
		Total	≤ 30mW		
Power dissipation can be reduced to 15 mW at 16 kHz sampling with 10-bit ADC resolution with a 20 MHz clk					
Digital I/Os	1.2V CMOS I/Os		Output impedance (fixed drive strength) ≤ 100 W		
	Outputs		Capacitive only loads		
			High-speed data / test	SPI output	
			≤ 10 pF, up to 160 MHz		≤ 40 pF, up to 40 MHz
	System inputs		rstb: active low reset input clk: system clock, 50% duty cycle, 160 MHz Rise-time = fall-time ≤ 25% of clk period, 1.56ns		
	4-wire SPI		csn: active low chip select, generated at falling sclck sdin: serial data input, generated at falling sclck sclck: serial clk (40 MHz, 1/4 <sup>th</sup> of clk frequency) (Rise-time = fall-time ≤ 6.25 ns) sdout: serial data output, sampled at falling sclck (Load for sdout ≤ 40 pF)		
	Serial data ports		160 Mbps at 160 MHz system clock Load for data ports ≤ 10 pF		
			Port-0	data_ser_out<0> wclk_ser_out<0>	
			Port-1	data_ser_out<1> wclk_ser_out<1>	
			Port-2	data_ser_out<2> wclk_ser_out<2>	
	Port-3	data_ser_out<3> wclk_ser_out<3>			
Package	Flip-chip CSP with Cu bumps		4 mm x 4 mm		
	Bump array / pitch		28 x 26	724 Bumps (none on corners)	140 μm
	Cu bump composition		40 μm Cu + 20 μm SnAg		
	Die thickness		12 mil (~305 μm)		

Active electrodes	Active	Reference	Test	
	512	8	8	
Input capacitance	6.8 pF			
Panel count	4 panels			
	One panel has 128 active electrodes, 2 reference, and 2 test electrodes			
Neural interface elements (nixels)	Selectable input configuration			
	Differential or single-ended operation			
	LNA + S/H + Comparator			
LNA parameters	Parameter name	Min	Max	
	Gain, input referred noise	13 V/V, 27 $\mu$ V rms	200 V/V, 4.5 $\mu$ V rms	
	High-pass corner	3.8 Hz	460 Hz	
	Low-pass corner	1.45 kHz	14 kHz	
Sample-and-hold (S/H)	Ping-pong architecture, with dual analog memory			
Comparator	In-nixel digitization with comparator Analog front-end of single-slope ADC			
ADC parameters	Single-slope-ADC, distributed architecture			
	User selectable resolution, 10–16-bit in resolution, $\leq$ 32 kHz in sampling			
	ADC Mode-0: <i>Low resolution, high sample rate</i>	10-bit @ 32 kHz sampling, 40 MHz clk		
	ADC Mode-1: <i>High resolution, mid sample rate</i>	12-bit @ 16 kHz sampling, 80 MHz clk		
	ADC Mode-2: <i>High resolution, high sample rate</i>	12-bit @ 32 kHz sampling, 160 MHz clk		
	ADC Mode-3: <i>Very high resolution low sample rate</i>	14-bit @ 8 kHz sampling, 160 MHz clk		
	ADC Mode-4: <i>Ultra high resolution very low sample rate</i>	16-bit @ 2kHz sampling, 160 MHz clk		
Active electrodes	Configurations	Total	Spike	LFP
	Differential, all for spike	512	480	32
	Diff. spike, single-ended LFP	496	480	16
	Single-ended spike, single-ended LFP	256	240	16
Reference electrodes	2 external reference electrodes / panel	Spike and LFP nixels have separate external reference electrode connections		
	2 external reference electrodes / chip			
	1 external reference electrodes / panel	Single external reference electrode		
	Programmable internal reference	Using internal DACs as reference		
Analog drive	On-chip bias generation	Band-gap based programmable bias generation using voltage and current mode DACs		
		12-bit v-DACs (references and test signals) 7-bit i-DACs		
	Ramp generation for ADC	Programmable ramp generation for ADC with variable slope and reset parameters		
Digital drive	Flexible and programmable operation of the chip	Programmable static circuit configuration of analog and digital circuits in terms of connectivity		
		Programmable timing generation for nixels and ADCs to adjust resolution and conversion time		
		Command based operation of the chip		
		Address based programming over SPI		

Programming interface	Serial	4-wire SPI
		32-bit SPI Words
		8-bit command, 8-bit address, 16-bit data
		"48 sclk cycles" per SPI operation
Test outputs	Analog	2 Analog voltage test output pads to monitor on-chip generated biases and reference voltages (pseudo-differentially)
		Voltage test outputs can be used to overwrite internally generated biases and references, including the global routing resources for reference and test electrodes used in the nixel array
	Digital	1 test current output to monitor on-chip generated reference and i-DAC output current 3-bit digital test outputs to monitor internal digital signals can also be configured to work as general-purpose output pins controlled over SPI

## Electrical Interface

The NYX1-512 chip runs on dual supply voltages of 1.2 V for digital blocks and 2.5 V for analog blocks. It uses 16-bit digital data paths internally and has four serial data ports with a 16:1 serialization ratio to send out the digital recording results. Each port has a dedicated word clock to indicate the word boundaries in the serial data stream.

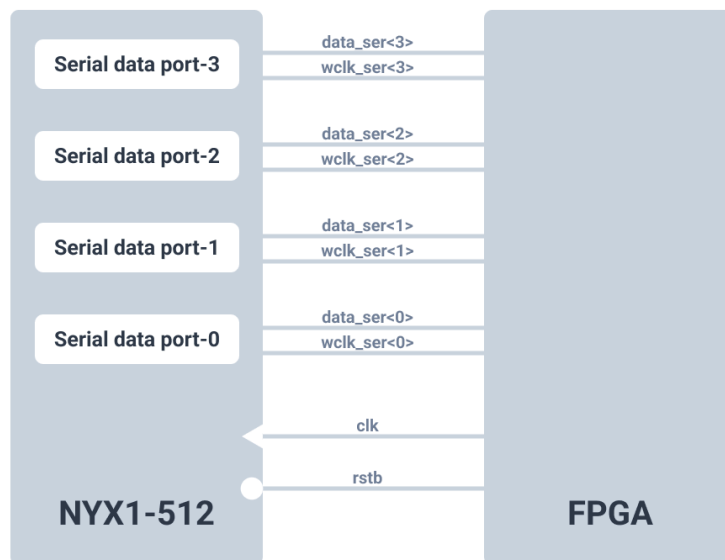


The NYX1-512 chip has five digital inputs and twelve digital outputs (I/Os) connected to the world using 1.2 V CMOS I/O buffers and powered with the 1.2 V digital supply of the chip. The chip also has 3-bit general purpose data outputs (GPOs) that can be controlled over SPI if needed.

### Digital I/Os


No.	I/O Name	Type	Specification
1	System	1.2V CMOS Input	Active low asynchronous reset
2			System clock, $\leq 160$ MHz, 50% duty cycle
3	4-wire SPI	1.2V CMOS Input	Active low chip select input for SPI
4			Serial data input for SPI
5			Serial clock for SPI, sclk freq = $\frac{1}{4}$ of clk
6			Serial data output of SPI
7-10	Serial data ports	1.2V CMOS Output	Serial data port for digitized nixels
11-14			Word clock for serial data port
15-17	Test / GPO	1.2V CMOS Output	Digital test outputs, can be hooked to SPI

The NYX1-512 chip uses a 4-wire SPI interface to program and control the chip, with active low chip select (csn), serial data input (sdin), serial clock (sclk), and serial data output (sdout). SPI interface uses 32-bit words, composed of an 8-bit command (<31:24>), an 8-bit address (<23:16>), and an 16-bit data (<15:0>) and sends the most-significant-bit (MSB) first. Since SPI operates at the rising edge of sclk, csn and sdin are applied at the falling edge of the sclk. Likewise, sdout from the SPI will be updated at the rising edge of the sclk, therefore it should be captured by the external electronics at the falling edge of sclk.

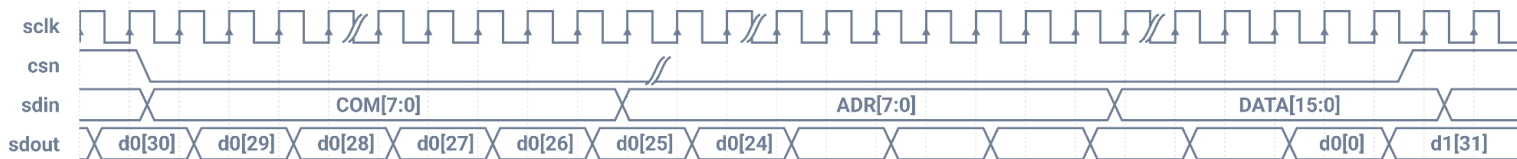


SPI timing takes 32 sclk cycles to enter the 32-bit SPI words into the input shift register of the SPI when csn is LOW. When csn is HIGH, it takes an additional 16 sclk cycles for the SPI controller to decode the SPI commands and read or write to the SPI registers. Including this idle time, an SPI operation will take 48 sclk cycles to complete. It is suggested to generate sclk from clk with a clock division ratio of 4.

Prior to streaming out a window of nixels, a panel keeps wclk and data low. For each enabled nixel in a panel (based on the window register settings), the panel will send a word. Each word is prefixed by a wclk high pulse—the corresponding 16 bits are sent out **starting when wclk goes low**—the data bits are valid on the *falling edge* of the clk:

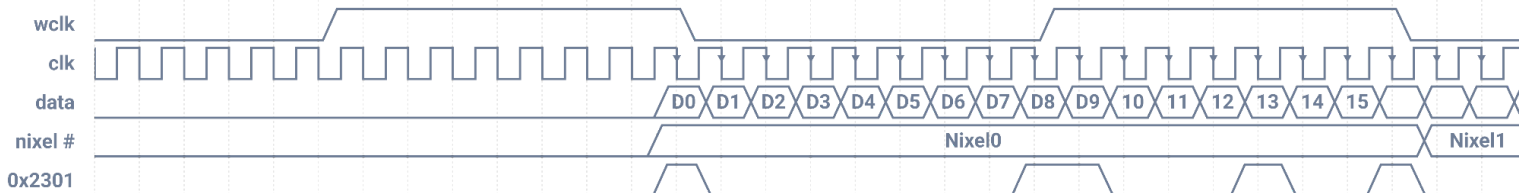

The last wclk will appear as:

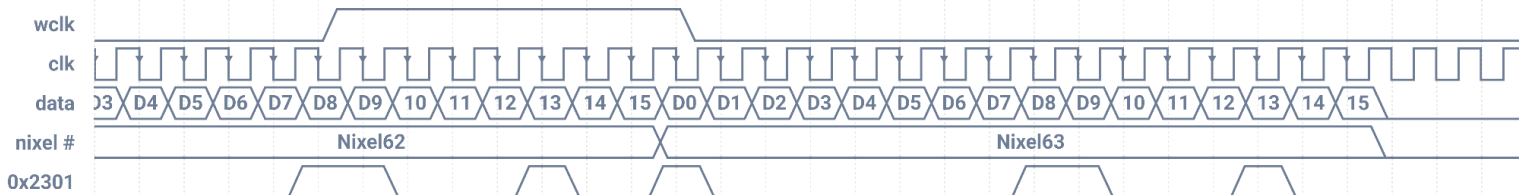

The NYX1-512 chip supports clk frequencies up to 160 MHz, providing ADC sampling rates up to 12-bit resolution at 32 kHz sampling rate. In this case, ADC sampling time will be 31.25  $\mu$ s, and the serial data stream for full resolution (64 nixels or 128 electrodes per panel) will only take 6.4  $\mu$ s, which is less than 21% of the available ADC sampling time. In cases where 10-bit ADC resolution will be enough, a 40 MHz clk will be sufficient for the same 32 kHz ADC sampling. In this 32 kHz 10-bit ADC mode, serial data can be transmitted in 25.6  $\mu$ s, which will correspond to about 82% of the available ADC sampling time. At clock frequencies less than ~33 MHz, it will not be possible to maintain 32 kHz ADC sampling rate without decreasing the ADC resolution below 10-bit. If lower resolution is not an option, then the ADC sampling rate can be reduced to 16 kHz or below to allow enough clock cycles for the single-slope type ADCs to perform ADC conversion. For example, 16 kHz ADC sampling with 10-bit resolution will be possible with clock frequencies just above 17 MHz.

## Command Structure

All NYX1-512 SPI commands are 32-bit values composed of an 8-bit command or operation code (OpCode) <31:24>, an 8-bit register address <23:16>, and 16-bit data or payload <15:0>. Below are the available operation codes followed by examples of common register operations.

OpCode	Name	Description	Expected behavior
0x00	NOP	No Operation	None
0x01	SOFT_RST	Soft Reset	Resets all registers to default values Requires 13 ms delay after reset
0x02	SOFT_CMD	Special Operation	Executes special commands
0x40	READ	Read Register	Returns register value
0xC0	WRITE	Write Register	Writes value to register

### Read from the DAC register

```
C/C++
uint32_t read_cmd = 0x40000000;           // OpCode for READ
uint32_t reg_addr = 0x02;                // Register address to read
uint32_t read_reg_cmd = read_cmd | (reg_addr << 16);
// Result: 0x40020000
// Expected response: 0x40020xxx where xxx is the register value
```

Register read operations return 16-bit values and require waiting for a response. To validate the response, you can use any of the following methods:

- Upper 8 bits of response should match sent OpCode
- Register address in response should match sent address
- Behavior is undefined for invalid register addresses

### Write to the DAC register

```
C/C++
uint32_t write_cmd = 0xC0000000;         // OpCode for WRITE
uint32_t dac_addr = 0x02;                // Register address (e.g., VDACC0_VCM_LNA)
uint32_t dac_value = 0x0908;            // Value to write
uint32_t write_dac_cmd = write_cmd | (dac_addr << 16) | dac_value;
// Result: 0xC0020908
```


Register write operations must provide 16-bit values and do not generate a response. DAC writes common use values between 0x0000 and 0x0FFF (12-bit DAC) while IDAC writes common use values between 0x00 and 0x7F (7-bit IDAC).

Both read and write operations require a valid register address to function properly (0x00 to 0x65).

## The Nixel Array

The nixel array is a 2D array of 256 neural recording elements, called nixels, which can each record from two electrodes simultaneously. Every nixel has an input switch matrix, a capacitively coupled fully-differential low-noise amplifier (LNA), an analog sample and hold (S/H) circuit, and an analog comparator.

The analog drive generates the required currents and voltages, as well as the static and dynamic reference and test signals used by the Nixel 512 chip. The digital drive generates all the control and timing signals used to configure and operate the chip using a 4-wire serial programming interface (SPI).

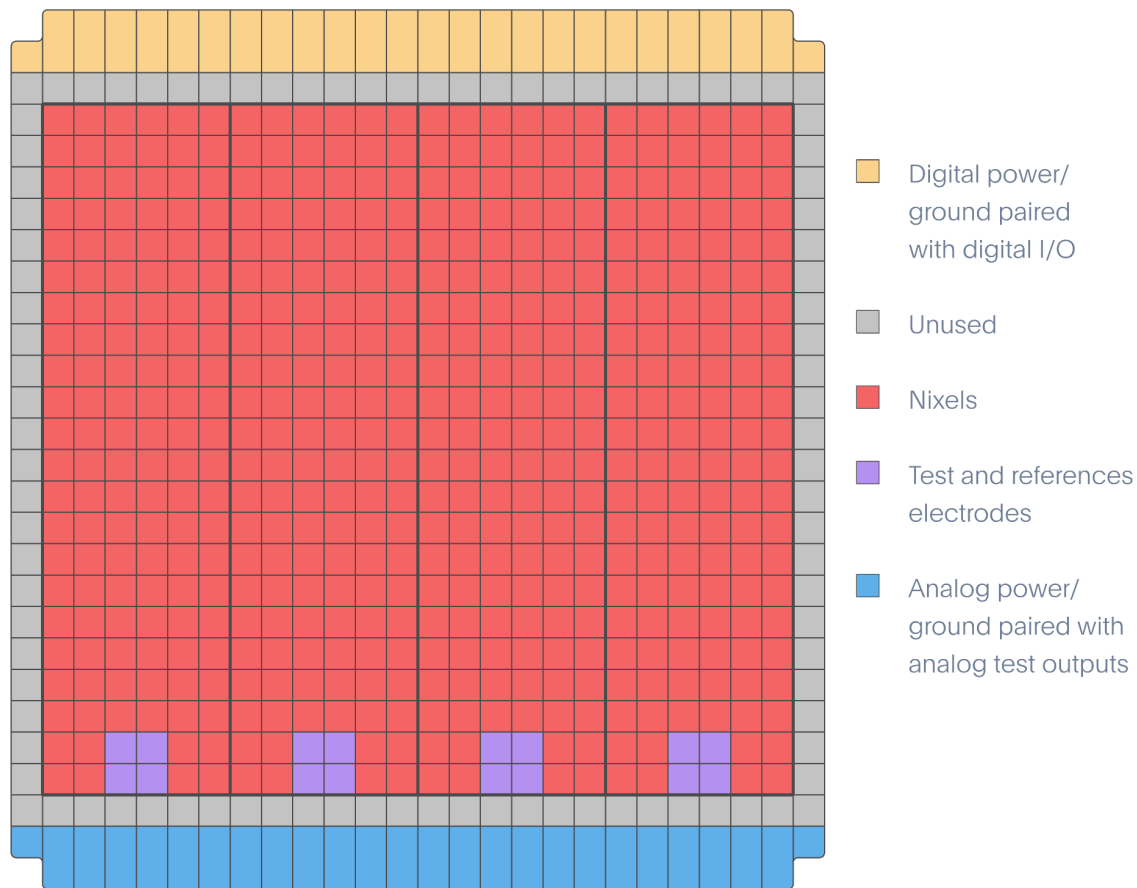

### Electrode Input Arrangement

The nixels are arranged in four identical panels, with each panel containing 128 electrode inputs. Each nixel is a differential amplifier connected to two electrode sites, labeled as even and odd.

The bottom two rows of nixels contain reference and test electrode inputs. Pin R1 is the LFP nixel reference, pin R2 is the spike nixel reference, and pins T1 and T2 can be used for testing purposes. The four adjacent nixels [0–7] are used for LFP recording while nixels [8–127] are used for spike recording. You can program the LFP nixels and spike nixels to have independent input switch configurations, gain, high-pass and low-pass filter corner settings.

All four panels can operate in parallel to record from all 512 electrodes at once. For the simplest configuration, you can connect all 16 reference and test electrode inputs together to a large reference electrode located in your active area.

122	123	124	125	126	127
116	117	118	119	120	121
110	111	112	113	114	115
104	105	106	107	108	109
98	99	100	101	102	103
92	93	94	95	96	97
86	87	88	89	90	91
80	81	82	83	84	85
74	75	76	77	78	79
68	69	70	71	72	73
62	63	64	65	66	67
56	57	58	59	60	61
50	51	52	53	54	55
44	45	46	47	48	49
38	39	40	41	42	43
32	33	34	35	36	37
26	27	28	29	30	31
20	21	22	23	24	25
14	15	16	17	18	19
8	9	10	11	12	13
4	5	T1	T2	6	7
0	1	R1	R2	2	3

## Recording Parameters

Panel Organization	Mode	Resolution		
	Panel 0	Nixels 0–63		
	Panel 1	Nixels 64–127		
	Panel 2	Nixels 128–191		
	Panel 3	Nixels 192–255		
ADC Configuration	Mode	Resolution	Sampling Rate	Minimum Clock
	0	10-bit	32 kHz	40 MHz
	1	12-bit	16 kHz	80 MHz
	2	12-bit	32 kHz	160 MHz
	3	14-bit	8 kHz	160 MHz
	4	16-bit	2 kHz	160 MHz
Bandwidth and Gain	High-pass corner	3.8 Hz to 460 Hz (s_Ina registers)		
	Low-pass corner	1.45 kHz to 14 kHz (s_Ina registers)		
	Gain settings	13 V/V (27 $\mu$ V rms noise) to 200 V/V (4.5 $\mu$ V rms noise)		

Each nixel corresponds to two electrodes: one odd, one even. The first four nixels of each panel are LFP nixels and the rest are spike nixels. If you need 512-channel operation, all nixels must be enabled.

### Troubleshooting

If you are not getting data:

1. Ensure the SPI clock is precisely 1/4 the main clock frequency
2. Verify proper power supply voltages (1.2V digital, 2.5V analog)
3. Check reference electrode configurations per panel

NOTE: The NYX1-512 ADC architecture uses a dual sample and hold configuration per channel. Due to a capacitor mismatch, there is a systematic alternating DC offset between readouts (ping-pong artifact). For high fidelity recordings, it is important to remove this error using a digital filter. This can be done in several ways, but in all cases the filter should have high rejection at half the sample rate.

## Recording Configuration

### Recording Setup Sequence

The NYX1-512 chip supports simultaneous recording from up to 512 electrodes using its four identical panels. This section describes the sequence of operations required to configure and start recording with the NYX1-512 ASIC.

### Reset

Reset the NYX1-512 chip by issuing the reset command (bring chip pin rstb low). Wait a minimum of 13 ms for the reset to complete.

```
C/C++
reset_nixel512();
// wait 13 ms for reset to take effect
```

### Enable Channels

Configure power-down registers. Nixels are organized in groups of 16 nixels per register, with a total of 16 registers (256 nixels total). Each nixel within a register can be individually enabled or disabled. This is done by writing 0 (enable) or 1 (power down) to registers 0x23 to 0x32.


```
C/C++
// enable all nixels except for first and last:
write_register(0x23, 0x0001);
write_register(0x24, 0x0000);
...
write_register(0x31, 0x0000);
write_register(0x32, 0x8000);
```

### Configure LNA Inputs

Each nixel has positive (LNA+) and negative (LNA-) low-noise amplifier inputs. Each input is connected to a switch that can be connected to a number of inputs detailed in the table below. Register 0x15 controls the switches for LFP nixels, and register 0x20 controls switches for the spike nixels. Each is a 16-bit register. Bits [15:11] set the positive input switch (spe[4:0]) and bits [10:6] set the negative input switch.

spe[4:0]	LNA+
0b0_0000	high-z
0b0_0001	sub
0b0_0010	ela<1>
0b0_0100	ela<0>
0b0_1000	elr_glb
0b1_0000	elt_glb<1>

sne[4:0]	LNA-
0b0_0000	high-z
0b0_0001	sub
0b0_0010	ela<0>
0b0_0100	ela<1>
0b0_1000	elr_glb
0b1_0000	elt_glb<0>

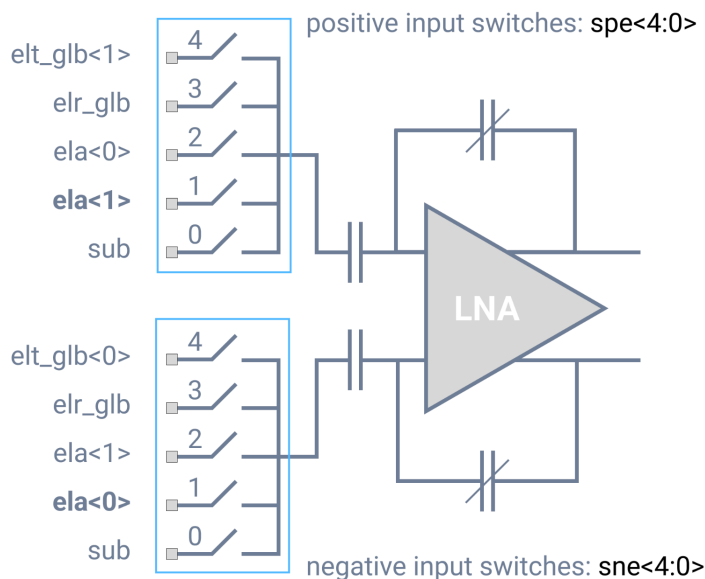

### Set Sample Rate and Bit Depth


The NYX1-512 chip uses a single slope ADC: a global counter and corresponding ramping signal both increment at the beginning of each sampling period. When the ramping signal surpasses the LNA's output voltage, the global counter is recorded and stored as that nixel's digitized, binary sample. The sampling rate is changed by controlling the stop value for the global counter (aka. line time) and ramping signal, as shown below:

```
C/C++
// Note: if using any of the on-chip DACs, we recommend using a line time that
// is a multiple of the DAC sampling rate
uint16_t new_line_time = (ASIC01_CLK_FREQ / desired_sampling_rate_u32) - 1;

// Ensure the same line time for all panels:
write_register(LINE_TIME_PANEL0, new_line_time);
write_register(LINE_TIME_PANEL1, new_line_time);
write_register(LINE_TIME_PANEL2, new_line_time);
write_register(LINE_TIME_PANEL3, new_line_time);
```

In the example above, bit depth is  $\log_2(\text{new\_line\_time})$ . E.g. for a 160 MHz clock and for 32 kS/s sampling rate, our line time is 5000 (for bit depth ~12).

In order to orchestrate proper timing of the ramp signal and counter, you must adhere to the following timing:

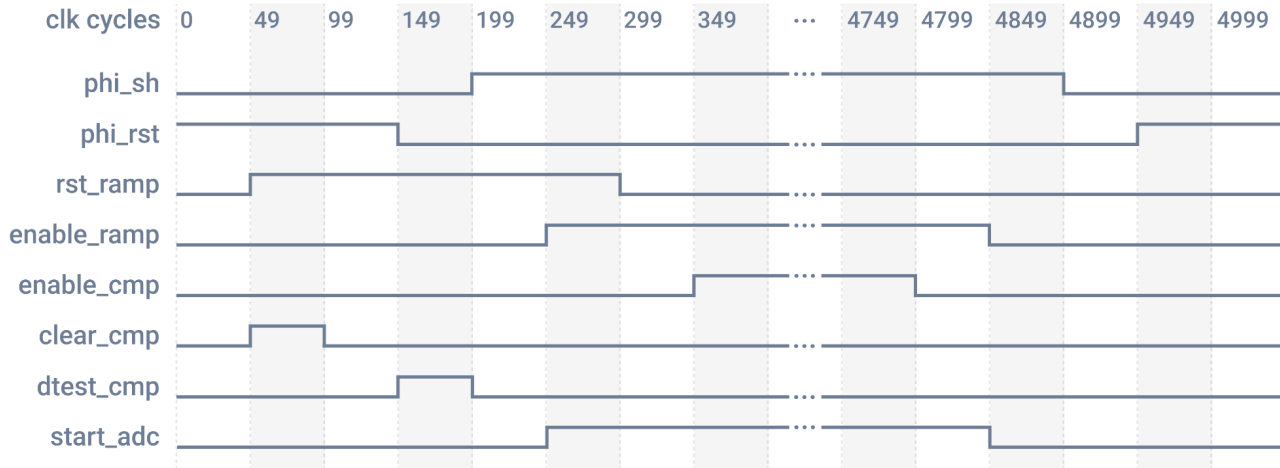

Internal Signals	SET	RESET
phi_sh	199	4899
phi_rstb*	149	4899
rst_ramp	49	299
enable_ramp	249	4849
enable_cmp	349	4799
clear_cmp	49	99
dtest_cmp	149	199
start_adc	249	4849
Line Time Duration	5000	

\*phi\_rstb - denotes the inversion of phi\_rst shown in the diagram above, where the reset is applied (the ramp is disabled) to start.

The example code below shows reasonable default values.

```
C/C++
// the global counter uses a 0-based index
write_register(RST_PHI_SH, new_line_time - 100);
write_register(RST_PHI_RSTB, new_line_time - 50);
write_register(RST_ENABLE_RAMP, new_line_time - 150);
write_register(RST_ENABLE_CMP, new_line_time - 200);
write_register(RST_START_ADC, new_line_time - 150);
```

In addition to the above global counter settings, you must also update the ramp signal appropriately. On chip DACs generate the ramp signal. You must ensure that the DAC starting and final voltage are in range.

Set the start voltage at 0.2 V or greater:

```
C/C++
double start_voltage = 0.5;
uint16_t new_ideal_start_code;
voltage_to_vdac_code(start_voltage, new_ideal_start_code);
write_register(VDAC04_VREF_INT, new_ideal_start_code);
```

Set the start voltage + ramp swing (`ramp_delta_v`) below 1.8V:

```
C/C++
uint16_t rst_enable_ramp = read_register(RST_ENABLE_RAMP);
uint16_t rst_rst_ramp = read_register(RST_RST_RAMP);
uint16_t integration_time_clocks = rst_enable_ramp - rst_rst_ramp;
double integration_time = integration_time_clocks /
    static_cast<double>(ASIC01_CLK_FREQ);
// integrator input current and integration cap set elsewhere --see docs
double ramp_delta_v = integration_time * integrator_input_current /
    integration_cap;
```

Set the ramp midpoint:

```
C/C++
double final_voltage = start_voltage + ramp_delta_v;
double ideal_midpoint_v = (start_voltage + final_voltage) / 2.0;
write_register(VDAC05_VREF_RAMP, new_ideal_midpoint_code);
```

Finally, set the sample rate to the following:

```
C/C++
double actual_sample_rate = ASIC01_CLK_FREQ / (new_line_time);
```

## Configure Reference Routing

NYX1-512 reference routing is configured via the `d_elref[15:0]` register (0x1C). Each panel (*k*) has its own reference routing, which can be set to either internal, or external mode (these refer to reference source). External reference mode is when the panel receives two reference inputs from the panel's external reference bus (`elr[1:0]`). Internal reference mode disconnects external inputs, even if they are present. When using internal reference, signal inputs are either left floating, or connected to internal test signal generators.

Reference is selected using a 4-bit binary signal applied per panel (`elref_k[3:0]`); 2-bits for external reference, and 2-bits for internal reference. This selection is made using 2 bits per mode. 0000, 1100, and 0011 are the most relevant options for reference selection.

- 0000 leaves reference inputs floating in a given panel (*k*:0, 1, 2, or 3).
- 1100 connects external references to the panel's global reference bus.
- 0011 connects the internal test signal generators, vDAC-7 and vDAC-8, to the panel's global reference bus (`elr_glb<1:0>`).

<code>elref_k[3:0]</code>	reference source
0b0000	high-z
0b0011	{ <code>vdac7_vtest</code> , <code>vdac8_vref_test</code> }
0b1100	<code>elr[1:0]</code>

## Initiate Start Sequence

Initiate the recording start sequence using one of the following SPI command sequences:

```
C/C++
// Normal recording:
const std::vector<uint32_t> START_CMDS = {0x00000000, 0xC0000010, 0x02000F00,
0xC036000F, 0xC05D00FF, 0x00000000, 0x00000000, 0x00000000};

// Fixed pattern output (by default, 0x2301):
const std::vector<uint32_t> START_DESER_TEST = {0x00000000, 0xC0000010,
0xC03AF924, 0x02000F00, 0xC05D00FF, 0x00000000, 0x00000000, 0x00000000};

// Incrementing output:
const std::vector<uint32_t> START_INCR_TEST = {0x00000000, 0xC0000010,
0xC03AF249, 0x02000F00, 0xC05D00FF, 0x00000000, 0x00000000, 0x00000000};
```

## Stop Recording

The sequence is stopped by issuing the reset command: `reset_nixel512()`;

Pin Specifications

NYX1-512 Chip			Center of Bump			I/O Type	Explanation	Value	Comment	Regulators	Direction	Avg. Current, Range, Resolution	
Rows	Col	Bump Name	X (um)	Y0 (um)	Y1 (um)								
A	B	1	sub	240	100	240	Analog power / ground	Substrate	0.0 V				
		2	vssa_bg	380				Ground / analog / bandgap	0.0 V				
		3	vdda_bg	520				Supply / analog / bandgap	2.5 V	1.8-2.8 V adjustable	REG-1	SOURCE	≤ 5 mA (DC)
		4	vdda_lna	660				Supply / analog / LNA	2.5 V		REG-2	SOURCE	≤ 10 mA (close to DC)
		5	vssa_lna	800				Ground / analog / LNA	0.0 V				
		6	vssa_cmp	940				Ground / analog / comparator	0.0 V				
		7	vdda_cmp	1080				Supply / analog / comparator	2.5 V		REG-3	SOURCE	≤ 10 mA (close to DC)
		8	dvdd2p5	1220				Supply / 2.5 V digital Blocks	2.5 V		REG-4	SOURCE	≤ 5 mA (switching)
		9	dvdd	1360				Supply / digital / core + IOs	1.2 V		REG-5	SOURCE	≤ 5 mA (switching)
		10	dvss	1500				Ground / digital / core + IOs	0.0 V				
		11	sub	1640				Substrate	0.0 V				
		12	vdda_lna	1780				Supply / analog / LNA	2.5 V				
		13	vssa_lna	1920				Ground / analog / LNA	0.0 V				
		14	vssa_cmp	2060				Ground / analog / comparator	0.0 V				
		15	vdda_cmp	2200				Supply / analog / comparator	2.5 V				
		16	dvdd2p5	2340				Supply / 2.5V digital blocks	2.5 V				
		17	dvdd	2480				Supply / digital / core + IOs	1.2 V				

NYX1-512 Chip			Center of Bump			I/O Type	Explanation	Value	Comment	Regulators	Direction	Avg. Current, Range, Resolution
Rows	Col	Bump Name	X (um)	Y0 (um)	Y1 (um)							
A	B	18	dvss	2620	100	240	Analog power / ground	Ground / digital / core + IOs	0.0 V			
		19	sub	2760				Substrate	0.0 V			
		20	vdda_lna	2900				Supply / analog / LNA	2.5 V			
		21	vssa_lna	3040				Ground / analog / LNA	0.0 V			
		22	vssa_cmp	3180				Ground / analog / comparator	0.0 V			
		23	vdda_cmp	3320				Supply / analog / comparator	2.5 V			
		24	itest_out	3460				Current test pad / analog / bias generator	0-32 μA sourcing (p-mirror)			
		25	vtest_out <0>	3600				Voltage test pad [0] / analog / bias generator	0-2.5 V v-DAC outputs			
		26	vtest_out <1>	3740				Voltage test pad [1] / analog / bias generator	0-2.5 V v-DAC outputs			
AG	AH	1	sub	240	3740	3880	Digital power / ground / IOs	Substrate	0.0 V			
		2	dvss	380				Ground / digital / core + IOs	0.0 V			
		3	dvdd	520				Supply / digital / core + IOs	1.2 V			
		4	csn	660				CMOS (1.2) input: SPI chip select bar	0-1.2 V	From FPGA, at negedge sclk, 50 Ω series res		
		5	sdin	800				CMOS (1.2) input: SPI data in	0-1.2V	From FPGA, at negedge sclk, 50 Ω series res		

NYX1-512 Chip			Center of Bump			I/O Type	Explanation	Value	Comment	Regulators	Direction	Avg. Current, Range, Resolution
Rows	Col	Bump Name	X (um)	Y0 (um)	Y1 (um)							
AG	AH	6	sclk	940	3740	3880	Digital power / ground / IOs	CMOS (1.2) input: SPI clock	0-1.2 V, fscclk = 1/4 fclk, ≤ 40 MHz, 50% DTC	From FPGA, at negedge sclk, 50 Ω series res		
		7	sdout	1080				CMOS (1.2) output: SPI data out	0 - 1.2 V	To FPGA, at posedge sclk, 50 Ω series res		
		8	rstb	1220				Active low asynch. reset	0 - 1.2 V	From FPGA, 50 or 20 Ω series res		
		9	clk	1360				Input: clock	0 - 1.2V, fclk ≤ 160 MHz, 50% DTC	From FPGA, 50 or 20 Ω series res		
		10	sub	1500				Substrate	0.0 V			
		11	dvss	1640				Ground / digital / core + IOs	0.0 V			
		12	dvdd	1780				Supply / digital / core + IOs	1.2 V			
		13	data_ser <0>	1920				CMOS (1.2) output: serializer data output [0]	SDR data rate, at rising edge clk, ≤ 160 Mbps	To FPGA, 50 or 20 Ω series res		
		14	wclk_ser <0>	2060				CMOS (1.2) output: serializer word clock output [0]	Frequency = 1/16 of SDR data rate	To FPGA, 50 or 20 Ω series res		

NYX1-512 Chip			Center of Bump			I/O Type	Explanation	Value	Comment	Regulators	Direction	Avg. Current, Range, Resolution
Rows	Col	Bump Name	X (um)	Y0 (um)	Y1 (um)							
AG	AH	15	data_ser <1>	2200	3740	3880	Digital power / ground / IOs	CMOS (1.2) output: serializer data output [1]	SDR data rate, at rising edge clk, ≤ 160 Mbps	To FPGA, 50 or 20 Ω series res		
		16	wclk_ser <1>	2340				CMOS (1.2) output: serializer word clock output [1]	Frequency = 1/16 of SDR data rate	To FPGA, 50 or 20 Ω series res		
		17	data_ser <2>	2480				CMOS (1.2) output: serializer data output [2]	SDR data rate, at rising edge clk, ≤ 160 Mbps	To FPGA, 50 or 20 Ω series res		
		18	wclk_ser <2>	2620				CMOS (1.2) output: serializer word clock output [2]	Frequency = 1/16 of SDR data rate	To FPGA, 50 or 20 Ω series res		
		19	data_ser <3>	2760				CMOS (1.2) output: serializer data output [3]	SDR data rate, at rising edge clk, ≤ 160 Mbps	To FPGA, 50 or 20 Ω series res		
		20	wclk_ser <3>	2900				CMOS (1.2) output: serializer word clock output [3]	Frequency = 1/16 of SDR data rate	To FPGA, 50 or 20 Ω series res		
		21	sub	3040				Substrate	0.0 V			
		22	dvss	3180				Ground / digital / core + IOs	0.0 V			
		23	dvdd	3320				Supply / digital / Core + IOs	1.2 V			

NYX1-512 Chip			Center of Bump			I/O Type	Explanation	Value	Comment	Regulators	Direction	Avg. Current, Range, Resolution
Rows	Col	Bump Name	X (um)	Y0 (um)	Y1 (um)							
AG	AH	24	digtest <0>	3460	3740	3880	Digital power / ground / IOs	CMOS (1.2) digital test output [0]	0-1.2 V	To FPGA, 50 or 20 Ω series res		
		25	digtest <1>	3600				CMOS (1.2) digital test output [1]	0-1.2 V	To FPGA, 50 or 20 Ω series res		
		26	digtest <2>	3740				CMOS (1.2) digital test output [2]	0-1.2 V	To FPGA, 50 or 20 Ω series res		

# On-Chip Registers

The NYX1-512 chip contains 101 writable (RAM) registers. Below are the descriptions and default values of all of the registers.

## AnaDrive

ADR	REG NAME	REG INFO / NET NAME	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
0	d_bgvref[7:0]	BandGap and Vref Settings	0000									test_bg[2:0]			start_bg	pd_bg	pd_ota_vref	pull_down_drv_vref	pull_up_drv_vref
1	d_irefdac[15:0]	Iref Settings	0840	pd_iref	d_iref[7:0]							d_idac[6:0]							
2	d_vdac0[12:0]	vdac0_vcm_lna	0908				pd_vdac_0	vdac_0[11:0]											
3	d_vdac1[12:0]	vdac1_vncas_lna	0908				pd_vdac_1	vdac_1[11:0]											
4	d_vdac2[12:0]	vdac2_vpcas_lna	0908				pd_vdac_2	vdac_2[11:0]											
5	d_vdac3[12:0]	vdac3_vcm_pre	0908				pd_vdac_3	vdac_2[11:0]											
6	d_vdac4[12:0]	vdac4_vref_int	073A				pd_vdac_4	vdac_4[11:0]											
7	d_vdac5[12:0]	vdac5_vref_ramp	0908				pd_vdac_5	vdac_5[11:0]											
8	d_vdac6[12:0]	vdac6_vimp	0172				pd_vdac_6	vdac_6[11:0]											
9	d_vdac7[12:0]	vdac7_vtest	0172				pd_vdac_7	vdac_7[11:0]											
10	d_vdac8[12:0]	vdac8_vref_test	0172				pd_vdac_8	vdac_8[11:0]											
11	d_vdac9[12:0]	vdac9_vncas	0908				pd_vdac_9	vdac_9[11:0]											
12	d_vdac10[12:0]	vdac10_vpcas	0908				pd_vdac_10	vdac_10[11:0]											
13	d_idac0[7:0]	idac0_vnbias_lna	0010									pd_idac_0	idac_0[6:0]						
14	d_idac1[7:0]	idac1_vpbias_pre	0010									pd_idac_1	idac_1[6:0]						
15	d_idac2[7:0]	idac2_vnbias_d2s	0010									pd_idac_2	idac_2[6:0]						
16	d_idac3[7:0]	idac3_i_int	0020									pd_idac_3	idac_3[6:0]						

AnaDrive

ADR	REG NAME	REG INFO / NET NAME	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0	
17	d_idac4[7:0]	idac4_ipbias_int	0020									pd_idac_4	idac_4[6:0]							
18	d_idac5[7:0]	idac5_ipbias_ramp	0020									pd_idac_5	idac_5[6:0]							
19	d_idac6[7:0]	idac6_inbias_vdac0	0050									pd_idac_6	idac_6[6:0]							
20	d_idac7[7:0]	idac7_inbias_vdac1_2	0050									pd_idac_7	idac_7[6:0]							
21	d_idac8[7:0]	idac8_inbias_vdac3	0050									pd_idac_8	idac_8[6:0]							
22	d_idac9[7:0]	idac9_inbias_vdac4	0050									pd_idac_9	idac_9[6:0]							
23	d_idac10[7:0]	idac10_inbias_vdac5	0050									pd_idac_10	idac_10[6:0]							
24	d_idac11[7:0]	idac11_inbias_vdac6_7_8	0050									pd_idac_11	idac_11[6:0]							
25	d_test[10:0]	Analog Test MUX Settings	0000							enable_i_test	sel_itest	enable_vtest	sel_vtest1[3:0]				sel_vtest0[3:0]			
26	d_ramp[5:0]	RampGen Settings	0006											ramp_test_enable	ramp_pd	ramp_scap[3:0]				
27	d_eltest[15:0]	Electrode MUX Settings: Test Electrodes	0000	eltest_3[3:0]				eltest_2[3:0]				eltest_1[3:0]				eltest_0[3:0]				
28	d_elref[15:0]	Electrode MUX Settings: Ref Electrodes	CCCC	elref_3[3:0]				elref_2[3:0]				elref_1[3:0]				elref_0[3:0]				
29	d_elbus[15:0]	Electrode MUX Settings: Test Read Busses	0000	elbus_3[3:0]				elbus_2[3:0]				elbus_1[3:0]				elbus_0[3:0]				
30	d_elimp[7:0]	Electrode MUX Settings: Impedance Read Busses	0000									elimp_3[1:0]		elimp_2[1:0]		elimp_1[1:0]		elimp_0[1:0]		

Nixel

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
31	s_sw[15:0]	LFP Input Switches	1080	spe_BOT[4:0]				sne_BOT[4:0]				simp_BOT[5:0]							
32	s_sw[31:16]	Spike Input Switches	1080	spe_TOP[4:0]				sne_TOP[4:0]				simp_TOP[5:0]							
33	s_lna[15:0]	LNA-BOT HP + LP Filter Corner Selection	0010	lp_BOT[3:0]			hp_BOT[7:0]							gain_cap_BOT[3:0]					
34	s_lna[31:16]	LNA-TOP HP + LP Filter Corner Selection	0010	lp_TOP[3:0]			hp_TOP[7:0]							gain_cap_TOP[3:0]					
35	pd[15:0]	Power-Down for 16 Nixels, Start Index = 0	FFFF	pd[15:0]															
36	pd[31:16]	Power-Down for 16 Nixels, Start Index = 16	FFFF	pd[31:16]															
37	pd[47:32]	Power-Down for 16 Nixels, Start Index = 32	FFFF	pd[47:32]															
38	pd[63:48]	Power-Down for 16 Nixels, Start Index = 48	FFFF	pd[63:48]															
39	pd[79:64]	Power-Down for 16 Nixels, Start Index = 64	FFFF	pd[79:64]															
40	pd[95:80]	Power-Down for 16 Nixels, Start Index = 80	FFFF	pd[95:80]															
41	pd[111:96]	Power-Down for 16 Nixels, Start Index = 96	FFFF	pd[111:96]															
42	pd[127:112]	Power-Down for 16 Nixels, Start Index = 112	FFFF	pd[127:112]															
43	pd[143:128]	Power-Down for 16 Nixels, Start Index = 128	FFFF	pd[143:128]															

Nixel

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
44	pd[159:144]	Power-Down for 16 Nixels, Start Index = 144	FFFF	pd[159:144]															
45	pd[175:160]	Power-Down for 16 Nixels, Start Index = 160	FFFF	pd[175:160]															
46	pd[191:176]	Power-Down for 16 Nixels, Start Index = 176	FFFF	pd[191:176]															
47	pd[207:192]	Power-Down for 16 Nixels, Start Index = 192	FFFF	pd[207:192]															
48	pd[223:208]	Power-Down for 16 Nixels, Start Index = 208	FFFF	pd[223:208]															
49	pd[239:224]	Power-Down for 16 Nixels, Start Index = 224	FFFF	pd[239:224]															
50	pd[255:240]	Power-Down for 16 Nixels, Start Index = 240	FFFF	pd[255:240]															
51	s_cmp[15:0]	Digital Output Mux Select for Nixel	AAAA	s_cmp_3_TOP[1:0]	s_cmp_3_BOT[1:0]	s_cmp_2_TOP[1:0]	s_cmp_2_BOT[1:0]	s_cmp_1_TOP[1:0]	s_cmp_1_BOT[1:0]	s_cmp_0_TOP[1:0]	s_cmp_0_BOT[1:0]								

ADC

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
52	N_start_adc[15:0]	Start Value for ADC Counter, glb setting	00FA	N_start_ADC[15:0]															
53	N_stop_adc[15:0]	Stop Value for ADC Counters, glb setting	1324	N_stop_ADC[15:0]															
54	{gray_adc[3:0], rstb_adc[3:0]}	Gray/Bin Count Modes + ADC_Reset_Bar	0000									gray_adc[3:0]				rstb_adc[3:0]			

DigCont

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0	
55	sel_panel[7:0]	Panel Select for Dig. Test Outs and Ramp Controls	0011									sel_panel_test_outputs[3:0]				sel_panel_ramp_control[3:0]				
56	sel_test[14:0]	Select Active Dig. Test Outputs, 4-bit x 3 Mux Select Signals, global for all Panels	7000	enable_digtest_out[2:0]			select_digtest_out_2[3:0]				select_digtest_out_1[3:0]				select_digtest_out_0[3:0]					
57	enable_oe_ser [11:0]	Enable Serializer for Panels, ON Signal for Data /Wclk for each Panel	0FFF				enable_serializer[3:0]				enable_wclk[3]	enable_data[3]	enable_wclk[2]	enable_data[2]	enable_wclk[1]	enable_data[1]	enable_wclk[0]	enable_data[0]		
58	enable_mode_select_hs[15:0 ]	Enable Horizontal_Scanner, Mode Select Horizontal_Scanner	F000	enable_hs[3:0]				test_mode_select_hs_3[3:0]		test_mode_select_hs_2[3:0 ]			test_mode_select_hs_1[3:0 ]			test_mode_select_hs_0[3:0 ]				
59	chip_port_id_hs[15:0]	8-bit Chip ID (default h00), and 4-Port IDs for Panels. 8-bit Port IDs defaults are: 11, 10, 01, 00 (default hE4)	00E4	chip_id[7:0]								port_id_3[1:0]		port_id_2[1:0]		port_id_1[1:0]		port_id_0[1:0]		
60	spi_scan_window_size[15:0]	Windowing Register for Panel-0: 8-bit Column Count, 8-bit Column Start	4000	column_count_0[7:0]								column_start_0[7:0]								
61	spi_scan_window_size[31:16 ]	Windowing Register for Panel-1: 8-bit Column Count, 8-bit Column Start	4000	column_count_1[7:0]								column_start_1[7:0]								
62	spi_scan_window_size[47:32 ]	Windowing Register for Panel-2: 8-bit Column Count, 8-bit Column Start	4000	column_count_2[7:0]								column_start_2[7:0]								
63	spi_scan_window_size[63:48 ]	Windowing Register for Panel-3: 8-bit Column Count, 8-bit Column Start	4000	column_count_3[7:0]								column_start_3[7:0]								
64	spi_scan_start_time[15:0]	Horizontal Scanner Start Time for Panel-0	0032	scan_start_time_0[15:0]																

DigCont

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
65	spi_scan_start_time[31:16]	Horizontal Scanner Start Time for Panel-1	0032	scan_start_time_1[15:0]															
66	spi_scan_start_time[47:32]	Horizontal Scanner Start Time for Panel-2	0032	scan_start_time_2[15:0]															
67	spi_scan_start_time[63:48]	Horizontal Scanner Start Time for Panel-3	0032	scan_start_time_3[15:0]															
68	spi_scan_data[15:0]	SPI Scan Data: Static Test Data from SPI (for the Last MUX in Data Path, Default h2301, which is Prod ID for ASIC01	2301	spi_scan_data[15:0]															
69	enable_adc_timing_limit[7:0]	Enable for ADC Timing Generators in DigCont and Enable Count Limit for ADC Counters in DigRecord	00FF									enable_adc_timing[3:0]				enable_adc_count_limit[3:0]			
70	dtest_mode_cmp[7:0]	8-bit Mode select for D-test Signal Generation for Nixel Comps. 3-bit per Panel.	00FF									dtest_mode_cmp_3[1:0]		dtest_mode_cmp_2[1:0]		dtest_mode_cmp_1[1:0]		dtest_mode_cmp_0[1:0]	
71	dtest_data_cmp[11:0]	12-bit dtest_data from SPI Register	0000					dtest_data_cmp_3[2:0]			dtest_data_cmp_2[2:0]			dtest_data_cmp_1[2:0]			dtest_data_cmp_0[2:0]		
72	spi_direct_control[15:0]	16-bit SPI Direct Control Register for all the on-chip generated timing signals, mainly for testing purposes	0000	enable_spi_direct_control[4:0]				start_adc_direct	enable_cmp_direct	clear_cmp_direct	dtest_cmp_direct[2:0]				phi_sh_direct[2:0]		enable_ramp_direct	rst_ramp_direct	
73	spi_set_rst_ramp[15:0]	Timing: SET_RST_RAMP	0032	spi_set_rst_ramp[15:0]															
74	spi_reset_rst_ramp[15:0]	Timing: RESET_RST_RAMP	012C	spi_reset_rst_ramp[15:0]															
75	spi_set_enable_ramp[15:0]	Timing: SET_ENABLE_RAMP	00FA	spi_set_enable_ramp[15:0]															

DigCont

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
76	spi_reset_enable_ramp[15:0]	Timing: RESET_ENABLE_RAMP	12F2	spi_reset_enable_ramp[15:0]															
77	spi_set_phi_sh[15:0]	Timing: SET_PHI_SH	00C8	spi_set_phi_sh[15:0]															
78	spi_reset_phi_sh[15:0]	Timing: RESET_PHI_SH	1324	spi_reset_phi_sh[15:0]															
79	spi_set_phi_rstb[15:0]	Timing: SET_PHI_RSTB	0096	spi_set_phi_rstb[15:0]															
80	spi_reset_phi_rstb[15:0]	Timing: RESET_PHI_RSTB	1356	spi_reset_phi_rstb[15:0]															
81	spi_set_dtest_cmp[15:0]	Timing: SET_DTEST_CMP	0096	spi_set_dtest_cmp[15:0]															
82	spi_reset_dtest_cmp[15:0]	Timing: RESET_DTEST_CMP	00C8	spi_reset_dtest_cmp[15:0]															
83	spi_set_clear_cmp[15:0]	Timing: SET_CLEAR_CMP	0032	spi_set_clear_cmp[15:0]															
84	spi_reset_clear_cmp[15:0]	Timing: RESET_CLEAR_CMP	0064	spi_reset_clear_cmp[15:0]															
85	spi_set_enable_cmp[15:0]	Timing: SET_ENABLE_CMP	015E	spi_set_enable_cmp[15:0]															
86	spi_reset_enable_cmp[15:0]	Timing: RESET_ENABLE_CMP	12C0	spi_reset_enable_cmp[15:0]															
87	spi_set_start_adc[15:0]	Timing: SET_START_ADC	00FA	spi_set_start_adc[15:0]															
88	spi_reset_start_adc[15:0]	Timing: RESET_START_ADC	12F2	spi_reset_start_adc[15:0]															
89	spi_line_time[15:0]	LINE TIME REGISTER for Panel-0. It is suggested to have same LINE time for all Panels.	1388	spi_line_time[15:0]															
90	spi_line_time[31:16]	LINE TIME REGISTER for Panel-1. It is suggested to have same LINE time for all Panels.	1388	spi_line_time[31:16]															

DigCont


ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
91	spi_line_time[47:32]	LINE TIME REGISTER for Panel-2. It is suggested to have same LINE time for all Panels.	1388	spi_line_time[47:32]															
92	spi_line_time[63:48]	LINE TIME REGISTER for Panel-3. It is suggested to have same LINE time for all Panels.	1388	spi_line_time[63:48]															
93	{enable_line_time[3:0], auto_line_time[3:0]}	8-bit ENABLE_TRIGGER_LINE_TIME_REGISTER: Enabled all Panales by Default, Expects SPI AUTO Trigger	00F0									enable_line_time[3:0]				auto_line_time[3:0]			
94	{enable_rstb_retime[4:0], enable_signal_retime[3:0]}	9-bit ENABLE_RSTB_SIGNAL_REGISTER. All RSTB and SIGNAL Retime functions are ON by Default.	01FF									enable_rstb_retime[4:0]				enable_signal_retime[3:0]			
95	spi_gpout[2:0]	3-bit SPI_GPOUT REGISTER. Use Reg#55 to configure to use SPI_GPOUT for digtest_out[2:0]	0000													spi_gpout[2:0]			


Unused

ADR	REG NAME	REG INFO	DEF (HEX)	15	14	13	12	11	10	9	8	7	6	5	4	3	2	1	0
96	unused_register0[15:0]	Can be used for SPI R/W Testing	0000	unused_register0[15:0]															
97	unused_register1[15:0]	Can be used for SPI R/W Testing	0000	unused_register1[15:0]															
98	unused_register2[15:0]	Can be used for SPI R/W Testing	0000	unused_register2[15:0]															
99	unused_register3[15:0]	Can be used for SPI R/W Testing	0000	unused_register3[15:0]															
100	unused_register4[15:0]	Can be used for SPI R/W Testing	0000	unused_register4[15:0]															
101	unused_register5[15:0]	Can be used for SPI R/W Testing	0000	unused_register5[15:0]															


## Packaging

The NYX1-512 chip measures 4 mm x 4 mm x 0.3 mm and uses chip scale packaging with Cu bumps. The 2D bump array counts 28 rows and 26 columns with 724 total bumps (no bumps on corners) and a uniform bump pitch of 140 μm.

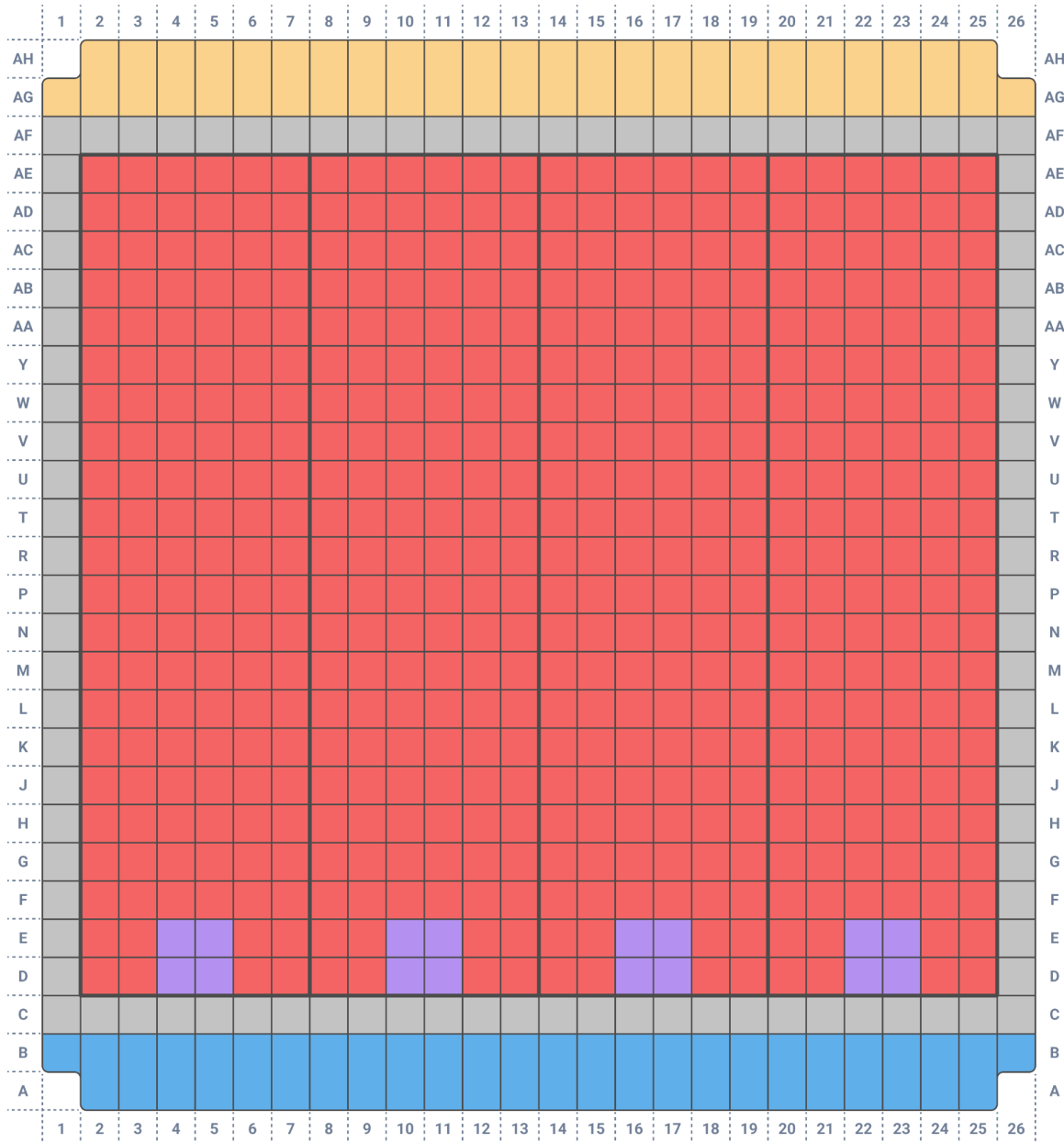

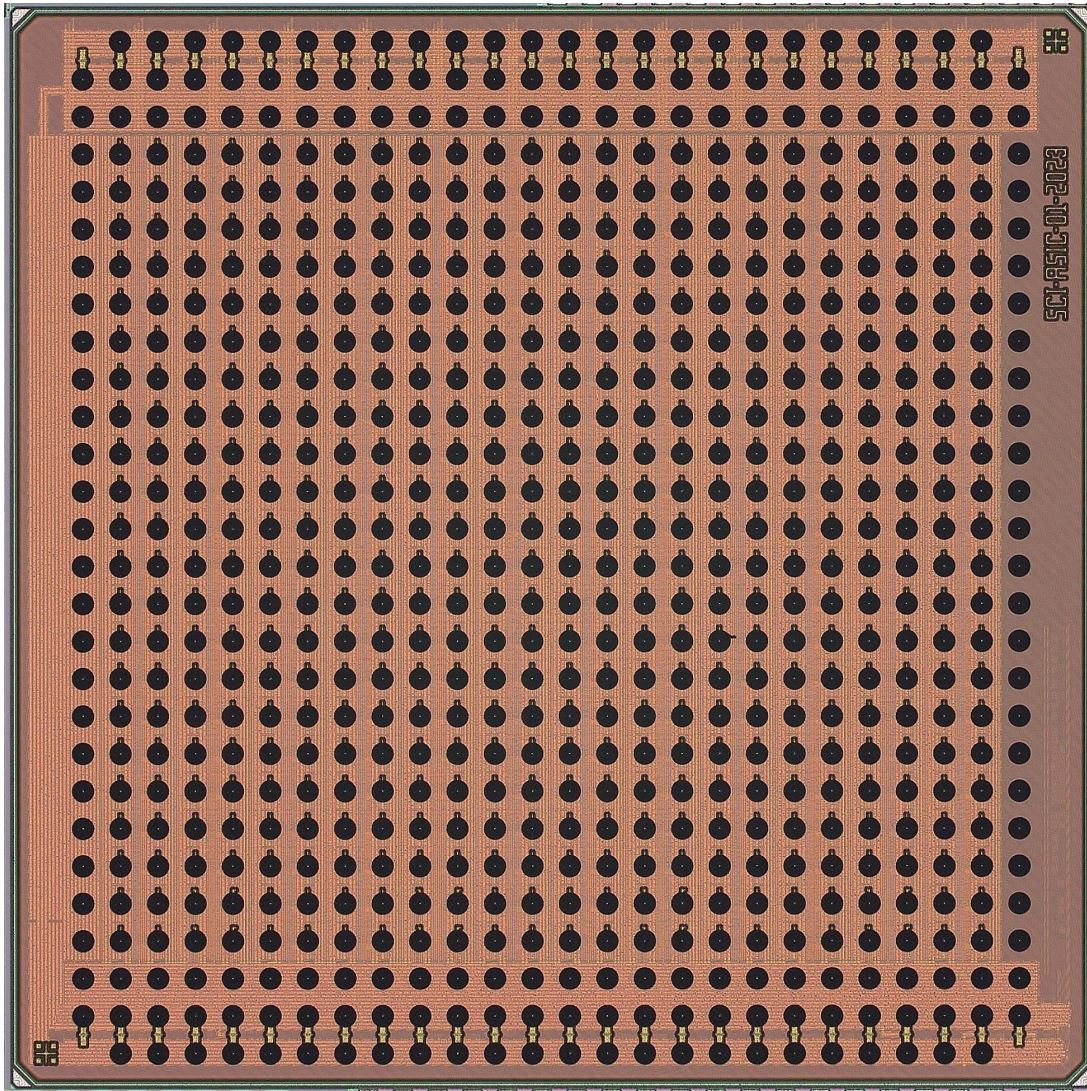

## Contact Information

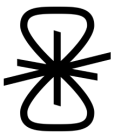

Science Corporation  
300 Wind River Way  
Alameda, CA 94501  
<https://science.xyz/support/>

*This preliminary product datasheet may change without notice.*